


Chunking Worms


1. Make sure the bottom of the plate to which you are moving the worms is labeled with the strain name and date.
2. Have both your old plate (starved worms, no bacteria) and new plate (no worms, bacteria) ready:

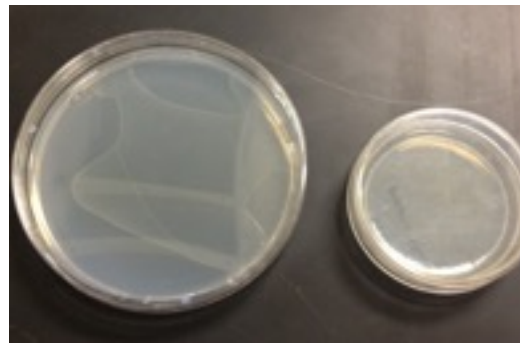

New Plate

Old Plate

3. Have the necessary tools ready:

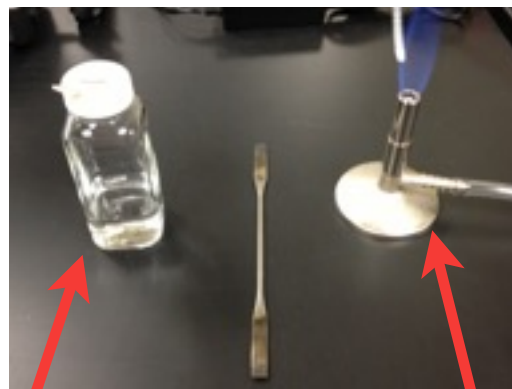


Jar of 95% Ethanol

Stainless Steel Spatula

Lit Bunsen Burner

4. Place the spatula in the jar of ethanol and then pass the spatula through the flame to sterilize it.

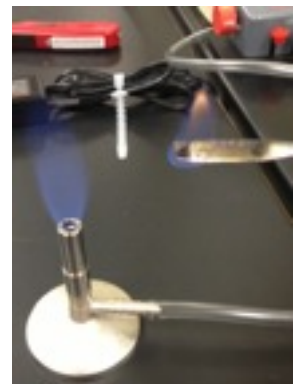

5. After the flame has extinguished from the spatula, use the sterilized spatula to cut through the agar in the center of the plate and cut out a $\sim 1 \text{ cm}^2$ area:



Making the first cut




A $\sim 1 \text{ cm}^2$ area

6. Remove the cut-out area using the spatula:



7. Place the agar chunk from the old plate **worm-side down** at the edge of the bacterial lawn on the new plate:

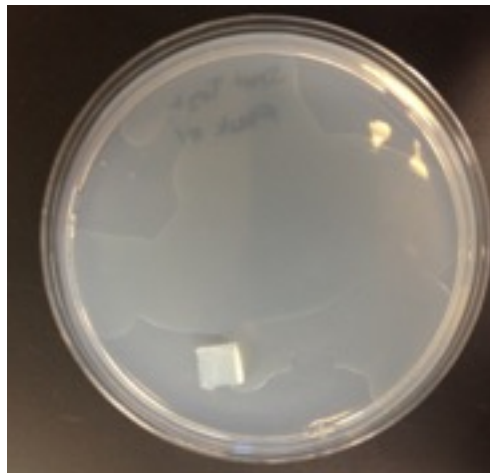

8. After you place the agar chunk down, hold the spatula in the flame of the Bunsen burner to melt/ burn any remaining agar and worms:



9. Place the spatula back in the ethanol to sterilize again before preparing the next sample.